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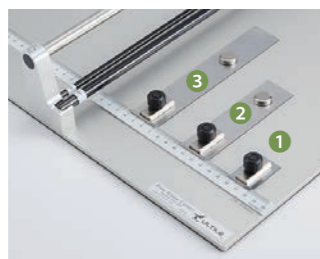
## 1. Magnetic substrate guide

### Standard | MG MGL MGU

Standard magnetic guides hold a substrate from both sides. Three sizes are available for various substrate sizes.


#### 1 MG 2 MGL 3 MGU

### Strong | MGA MGLA MGUA


Strong magnetic guides are used for cutting substrates repeatedly in the same size. Use these guides with the standard magnetic guides.
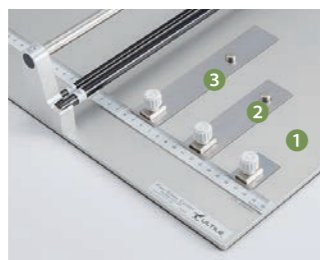
#### 1 MGA 2 MGLA 3 MGUA

### Precision | MGP

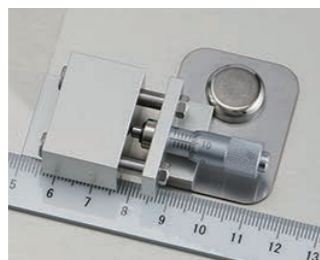The precision magnetic guide enables the user to cut in an accuracy less than 0.1mm. Using a caliper, it can further adjust and cut the substrate in an accuracy of  $\pm 30\mu\text{m}$ .

## 4. Option blade

### Carbide cutter

The carbide cutter is suitable for cutting standard hardness materials such as soda-lime glass and silicon substrates. ※-1. 2. 3

type	Normal			High penetrating
Blade	Carbide wheel for ultra-thin substrate	Carbide wheel for thin substrate	Carbide wheel for standard substrate	Carbide laser wheel
Thicknesses of target substrate	0.15 ~ 0.5mm	0.5 ~ 1.3mm	1.1 ~ 3.0mm	0.05 ~ 2.0mm
Color / stamp	Silver · U	Silver · T	Silver · N	Silver · LASER

### Diamond cutter

The diamond cutter has a long life and can cut a wide range of materials from standard hardness materials such as glass to very hard materials such as sapphire. ※-1. 2. 3

type	Normal			High penetrating
Blade	Diamond wheel for ultra-thin substrate	Diamond wheel for thin substrate	Diamond wheel for standard substrate	Diamond laser wheel
Thicknesses of target substrate	0.15 ~ 0.5mm	0.5 ~ 1.3mm	1.1 ~ 3.0mm	0.05 ~ 2.0mm
Color / stamp	Black · U	Black · T	Black · N	Black · LASER

※-1 The stamp can be found near the threaded screw hole for fixing the wheel.

※-2 The thicknesses of target substrate is for the case of soda-lime glass.

※-3 A diamond laser wheel is a high-performance wheel applicable to a wide range of substrate thicknesses, although it is not suitable for getting smooth fracture surfaces.

## 2. Positioning Magnifier

### Dedicated magnifier | LP

This is a dedicated 5x magnifier. It can be used either with a holder for scale magnifier or a magnifier holder for checking cutting position.

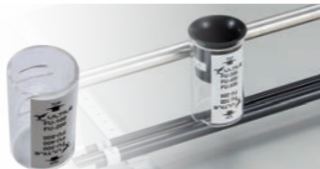
### Holder for scale magnifier | LHS

This holder enables the user to look at the scale perpendicularly. That makes substrate positioning easier.

### Magnifier holder for checking cutting position | LHC


This holder holds the magnifier to check the cutting position. Set the holder upside down when using FU-300 glass cutter or larger.

### Digital microscope system | DMS

Connect the microscope to your smartphone or computer to view the cut position on the screen at high magnification. Set the material on the attached microscope holder when using this system.

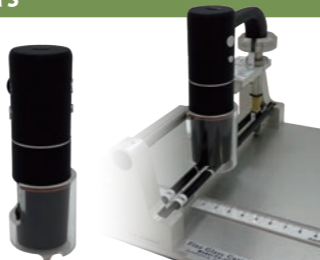※ Only available with FU-100 and FU-200.


## 3. Option parts for small-sized glass/substrate

### Slide holder | SSH-A

The slide holder is convenient for cutting substrates to very small sizes of 10 mm or less.

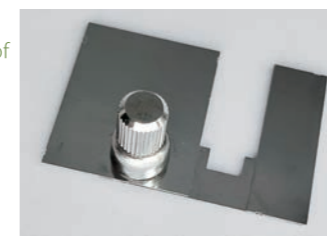※Compatible with substrates from 4 to 10mm.

Use magnetic substrate guides to hold the substrate from both ends.


The width of the right side portion of the slide holder is 10 mm. When cutting the substrate, place it in the slide holder and set the edge of the right slide holder at the target dimension +10mm.

※Example of cutting to 5mm

### Running tweezers | TSB

Just by pinching the substrate scribed by ULTILE, you can easily break micro-sized substrates less than 10mm while it is difficult to cut with your fingers. Furthermore, the substrate is held by the running tweezers even after it is cut, you don't have to worry about finding the top or bottom side of the substrate by dropping it.

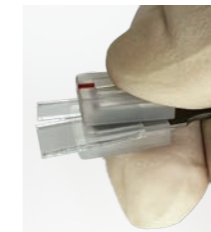We recommend this to break Si wafers.


## 5. Optional head

For cutting films

### Film cutting head | FCH

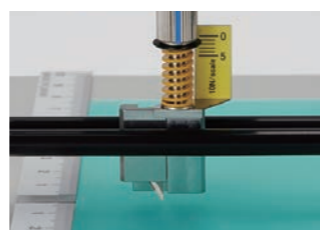This head can replace the wheel cutter. Use a dedicated mat for protecting the cutter from damage.

**Attaching a cutter** Set the cutter to the head. Put a cover for protecting the cutter from falling off.

### FCH dedicated cutter | FCC

Thirty spare cutters dedicated for the film cutting head.

### FCH dedicated mat | FCMS FCMM

A mat for cutting films Size S is the only mat usable with FU-100.

1 Size S 105×150 mm **FCMS**  
2 Size M 210×300 mm **FCMM**

For mechanical scriber

### Mechanical scriber head set | SBS

This head can replace the wheel cutter. A jig for adjusting the height of the scriber bit is included.

**Setting the bit** Place a sheet of thin paper at the bottom of the height adjustment jig where the bit touches it. Screw the bit and remove the head from the jig.

### Diamond bit | SBDA SBDB SBDC

Three sizes are available for various substrate sizes.

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
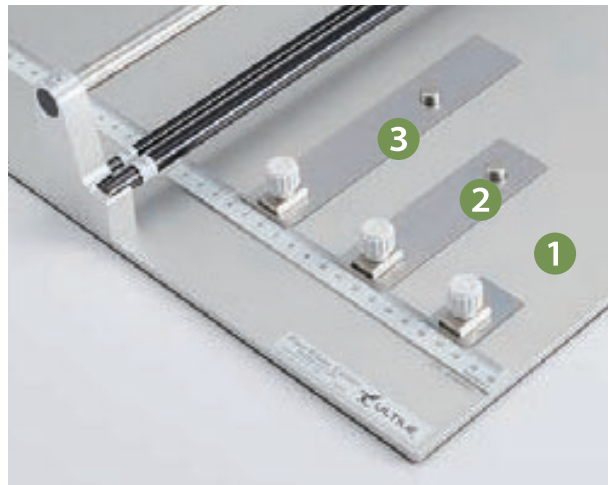
① MGA ② MGLA ③ MGUA

## Precision | MGP

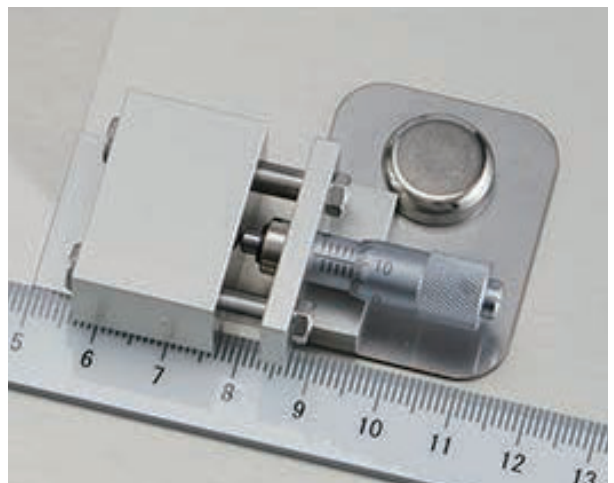The precision magnetic guide enables the user to cut in an accuracy less than 0.1mm. Using a caliper, it can further adjust and cut the substrate in an accuracy of  $\pm 30 \mu\text{m}$ .

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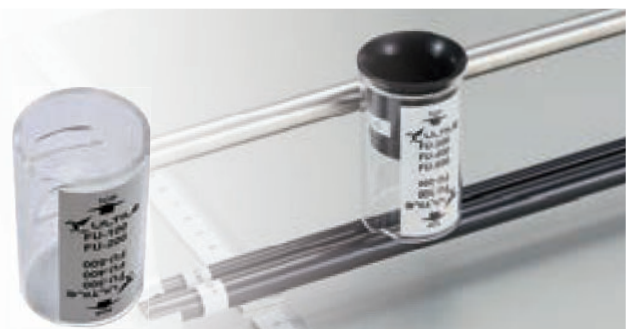
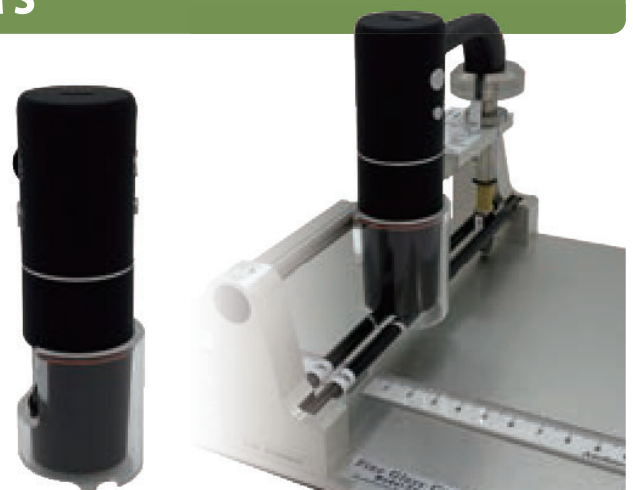
※ Only available with FU-100 and FU-200.



### 3.Option parts for small-sized glass/substrate

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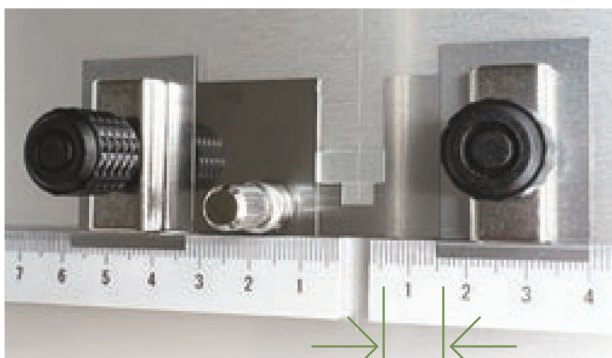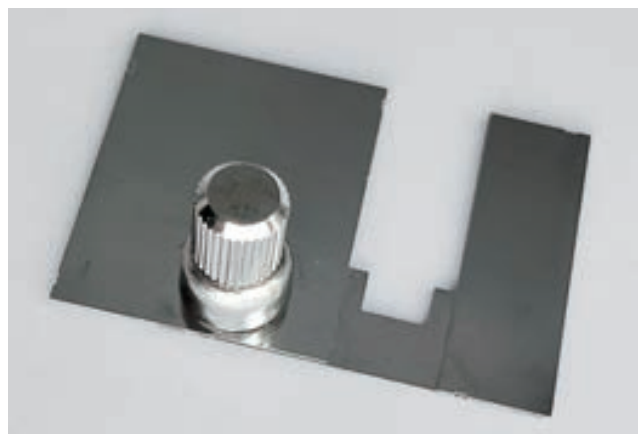## 4.Option blade

### Carbide cutter

The carbide cutter is suitable for cutting standard hardness materials such as soda-lime glass and silicon substrates. ※-1.2.3

type	Normal			High penetrating
Blade	Carbide wheel for ultra-thin substrate	Carbide wheel for thin substrate	Carbide wheel for standard substrate	Carbide laser wheel
Thicknesses of target substrate	0.15 ~ 0.5mm	0.5 ~ 1.3mm	1.1 ~ 3.0mm	0.05 ~ 2.0mm
Color / stamp	Silver • U	Silver • T	Silver • N	Silver • LASER

### Diamond cutter

The diamond cutter has a long life and can cut a wide range of materials from standard hardness materials such as glass to very hard materials such as sapphire. ※-1.2.3

type	Normal			High penetrating
Blade	Diamond wheel for ultra-thin substrate	Diamond wheel for thin substrate	Diamond wheel for standard substrate	Diamond laser wheel
Thicknesses of target substrate	0.15 ~ 0.5mm	0.5 ~ 1.3mm	1.1 ~ 3.0mm	0.05 ~ 2.0mm
Color / stamp	Black • U	Black • T	Black • N	Black • LASER

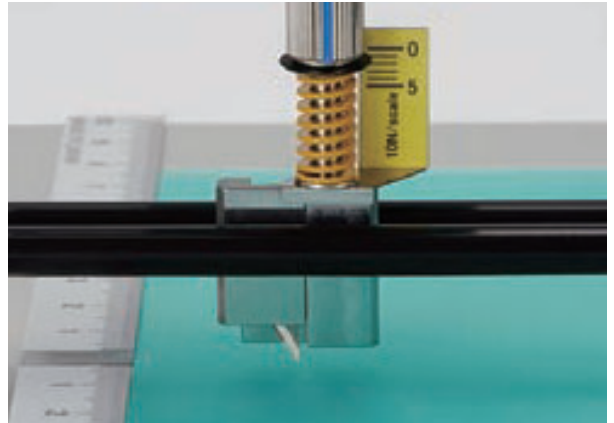
- ※-1 The stamp can be found near the threaded screw hole for fixing the wheel.
- ※-2 The thicknesses of target substrate is for the case of soda-lime glass .
- ※-3 A diamond laser wheel is a high-performance wheel applicable to a wide range of substrate thicknesses, although it is not suitable for getting smooth fracture surfaces.

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#### Attaching a cutter

Set the cutter to the head.

Put a cover for protecting the cutter from falling off.

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### FCH dedicated mat | FCMS FCMM

A mat for cutting films. Size S is the only mat usable with FU-100.

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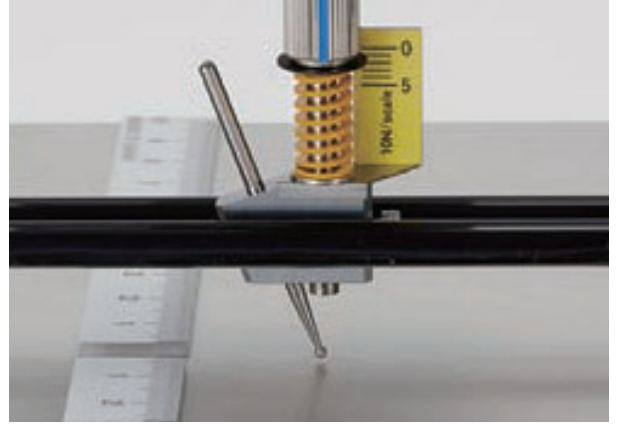
### Setting the bit

Place a sheet of thin paper at the bottom of the height adjustment jig where the bit touches it.

Screw the bit and remove the head from the jig.

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Astellatech, Inc.